



KNIFE MILL HM100

- Feed size □ <50mm
- Final fineness□□ 320μm

The Knife Mill HM100 is designed for processing soft, medium-hard, brittle, and fibrous samples, particularly those with high water, oil, or fat content.

APPLICATIONS

- Sample types: Soft, elastic, fibrous, water-containing, oily, fatty, dry samples.
- Examples of applications:
 - Agriculture: grains, feeds, etc.
 - Biology: mouse tissues, plant leaves, seeds, seedlings, and germs.
 - Medicine: traditional medicine, pills, etc.
 - Food: dried melons, vegetables, frozen food, candy, preserved fruits, and meat.

FEATURES

- Option for an autoclavable container
- Grinding container available in various materials
- Easy operation and quick startup
- Powerful 1.1KW motor with adjustable speed
- Achieve both coarse and fine grinding in a single step
- Three optional modes: interval mode, reverse mode, and press-and-play mode
- Equipped with overload protection and a safety lock design

THEORY & METHOD

Working Principle

The HM100 features two-blade knives positioned at the center of the grinding cup's bottom. These blades maintain a certain height gap between them. When powered by a 1,100W motor, the knives begin to rotate. The sharp steel knife edges cut effectively, ensuring reliable crushing and the creation of uniform phase effects. You can opt for the blunt side for coarse grinding of bulk samples or the tough side for fine grinding of samples.

The motor speed is electronically controlled and can be preset. This consistent speed ensures efficient and high-quality grinding homogenization. It takes just 10 to 30 seconds to process samples that are typically difficult to crush, yielding homogenized samples ready for analysis.

Knife Mill

TECHNICAL SPECIFICATIONS

Feed size	< 40mm
Final fineness	< 300μm
Speed	2000-10000rpm
Time setting	0 01-9 59 min/s
Batch quantity	150-700ml
Interval mode	Yes
Knives	stainless steel, titanium alloy

Number of blades	2
Rated power	1.1KW
Power supply	220V□ 50/60Hz (110V Also available upon request)
Instrument size	260*343*454mm
Package size	480*550*720mm
Net weight	16kg



Torontech Inc.

251 Consumers Road, Suite 1200
Toronto, Ontario, M2J 4R3, Canada 

Phone: +1 416 368 2721 | Fax: +1 416 981 7652
Email: info@torontech.com

Torontech, LLC

601 Brickell Key Drive, Suite 700
Miami, FL 33131, USA 

Toll-Free: 1-866-383-7919
Email: sales@torontech.com

Torontech ME FZE

Meydan Grandstand, 6th Floor,
Meydan Road, Nad Al Sheba,
Dubai-261440, UAE 

Phone: +971 488 19252
Email: me@torontech.com

