



GOLD PURITY DENSITY TESTER TTD-LK

Gold Purity Density Tester TTD-LK designed specifically for precious metals, accommodating general-size samples with dimensions of 210 x 170 x 120mm.

Maximum Weight Capacity: 1500g / 2000g / 3000g / 6000g

Weighing Precision: 0.01g / 0.01g / 0.01g / 0.1g

Purity Precision: 100% or 100.0%

Density Meter, Density Tester, Densimeter, Specific Gravity Tester

Density Meter, Density Tester, Densimeter, Specific Gravity Tester

Density Meter, Density Tester, Densimeter, Specific Gravity Tester

Density Meter, Density Tester, Densimeter, Specific Gravity Tester

APPLICATIONS

Suitable for

Ideal for use in the jewelry industry, banks, pawnshops, and gold and precious metal research laboratories.

Principle

Using the Archimedean principle's buoyancy method, gold karat, density, and purity percentage can be rapidly and accurately displayed.

Mode

1. Gold Purity Measurement Mode – Measures gold purity.
2. Platinum Purity Measurement Mode – Measures platinum purity.
3. Alloy Purity and Density Measurement Mode 1 – Measures the purity of other precious metals.
4. Alloy Purity and Density Measurement Mode 2 – Measures the purity of other precious metals.

FEATURES


- Gold Purity Density Tester TTD-LK equipped with RS-232 and USB interfaces for connectivity.
- Infrared temperature sensors automatically detect and compensate for water temperature variations.
- Enables quick and accurate determination of precious metal purity percentage and gold Karat number in just two measurement steps.
- Leaves no trace of contamination after the measurement, eliminating the need for sulfuric acid solution or standard test samples.
- Special software design directly displays gold Karat number and purity percentage, including variations like gold-copper-silver, gold-silver, and gold-copper.
- Features a setting function for the lower limit of the Karat value, allowing for Karat number adjustment from 1 to 24K.
- Offers a function to set the lower limit of PT value, allowing PT value adjustment from 1 to 1000.

- Capable of measuring the purity of various other precious metals, including platinum, silver, palladium, copper, rhodium, and more.
- Unable to test samples with gems and hollows; however, it can function as a general balance.
- Utilizes a large tank design to minimize errors due to the buoyancy of the hanging rail. The water tank size is 210 × 170 × 120mm.

TECHNICAL SPECIFICATIONS

Model		TTD-LK1500	TTD-LK2000	TTD-LK3000	TTD-LK6000
Max Weight		1500g	2000g	3000g	6000g
Weighing Precision		0.01g			0.1g
Density Precision		0.001g/cm ³			
Gold Mode	Karat	1K and 0.1K display (switchable)			
	Purity	1% and 0.1% display (switchable)			
PT Mode	PT	25PT and 1PT display (switchable)			
	Purity	1% and 0.1% display (switchable)			
Alloy Mode		1% and 0.1% display (switchable)			

Density Meter, Density Tester, Densimeter, Specific Gravity Tester

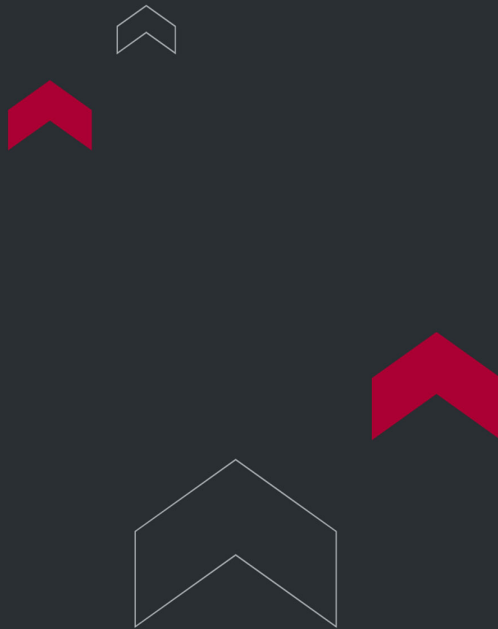

Torontech Inc.

251 Consumers Road, Suite 1200
Toronto, Ontario, M2J 4R3, Canada 

Phone: +1 416 368 2721 | Fax: +1 416 981 7652
Email: info@torontech.com

Torontech, LLC

601 Brickell Key Drive, Suite 700
Miami, FL 33131, USA 

Toll-Free: 1-866-383-7919
Email: sales@torontech.com

Torontech ME FZE

Meydan Grandstand, 6th Floor,
Meydan Road, Nad Al Sheba,
Dubai-261440, UAE 

Phone: +971 488 19252
Email: me@torontech.com

