



WET-RESISTANT BACTERIAL PENETRATION TESTER

Wet-resistant Bacterial Penetration Tester ToronFM™ WRBPT designed to evaluate how effectively a material resists bacterial penetration in liquids under mechanical friction. This device assesses the material's ability to block bacterial intrusion during frictional contact in a controlled chamber. Common applications include testing surgical gowns, medical apparel, and cleanroom garments

APPLICATIONS

Standards

1. ISO 22610-2018

Principle

The specimen was placed on an agar plate, followed by a fungus plate of the same size positioned on top. A high-density polyethylene film, approximately 10 microns thick, was then placed over the fungus plate. These three layers (with the test material at the bottom, the fungus plate in the middle, and the polyethylene film on top) were secured using a tapered steel ring.

The test involved subjecting the material to controlled movement across the surface of the petri dish, exposing it to combined pressure and friction. This simulation replicates real-world conditions where barrier materials face microbial penetration under wet environments.

During testing, microbes from the fungal plate penetrated the specimen and migrated to the surface of the agar medium. Colony counting was performed to quantitatively evaluate the material's resistance to microbial penetration.

TECHNICAL SPECIFICATIONS

Model	ToronFM™ WRBPT
Rotation Time of Rotating Table	Automatically controlled by the timer
Test Finger	Guided by the rotating outer wheel, can do lateral reciprocate operations from the center of the rotating
Test Finger Working Pressure to the Material	Adjustable
Material of the Test Parts	Corrosion-resistant stainless steel
Rotary Speed	(60 ± 1) RPM
Test Finger Pressure on the Material	(3 ± 0.02) N
The Rotation Speed of the Outer Wheel	5 ~ 6 RPM
The Timer Setting Range	0 ~ 99.99 min
Total Weight of Inner and Outer Ring	(800 ± 1) g



Torontech Inc.

251 Consumers Road, Suite 1200
Toronto, Ontario, M2J 4R3, Canada 

Phone: +1 416 368 2721 | Fax: +1 416 981 7652
Email: info@torontech.com

Torontech, LLC

601 Brickell Key Drive, Suite 700
Miami, FL 33131, USA 

Toll-Free: 1-866-383-7919
Email: sales@torontech.com

Torontech ME FZE

Meydan Grandstand, 6th Floor,
Meydan Road, Nad Al Sheba,
Dubai-261440, UAE 

Phone: +971 488 19252
Email: me@torontech.com

