



DESK SPUTTER & CARBON COATER - DSCR

The Desk Sputter & Carbon Coater - DSCR is a compact coating system that combines both sputtering and carbon coating capabilities. This versatile system can sputter noble metals like gold (Au), palladium (Pd), platinum (Pt), and gold/palladium (Au/Pd) onto non-conductive or poorly conductive specimens, and it can also apply carbon films to the samples.

APPLICATIONS

The Desk Sputter Coater model DSCR is specifically designed as a sample preparation system for Scanning Electron Microscopy (SEM). It is equipped with a rotary pump that enables the achievement of a vacuum level below 50 mTorr. This vacuum range is well-suited for sputter coating with noble metals and carbon fiber evaporation coating.

For sputtering deposition involving oxidizing metals, Field Emission SEM (FE-SEM), SEM with Energy Dispersive X-ray Spectroscopy (SEM EDX), and applications related to Transmission Electron Microscopy (TEM), we recommend considering other models such as DST1, DSCT, and DCT, which are better suited for these specific purposes.

Desk Sputter Coater

FEATURES

- Easy to operate
- Intuitive touchscreen for seamless control of the coating process and quick data input
- User-friendly software enabling both full-automatic and semi-automatic control
- Effortless loading and unloading of specimens
- Consistent film thickness depositions for repeatability
- Precise and reproducible coatings
- Capable of achieving high-resolution fine coatings (on the order of 2nm for Gold grain)

- Low voltage sputtering
- GLAD (Glancing Angle Deposition) Sputtering
- Plasma Cleaning
- No cooling system required
- Carbon coating available in both flash and pulse modes
- Dual carbon fiber for thicker layer deposition
- Covered by a two-year warranty for peace of mind
- CE conformity ensures compliance with safety standards
- Worldwide coverage for public and product liability insurance, offering protection in case of property damage or personal injury caused by Vac Coat systems.

Pulsed Carbon Fiber Evaporation

The small chamber of the desk vacuum sputter and carbon coater has the capability to perform pulsed carbon fiber evaporation. This method involves using short pulses for more precise and controlled deposition, effectively reducing the amount of debris typically associated with traditional carbon deposition processes.

Carbon Coating Head

The DSCR is equipped with an interchangeable head designed for carbon coating purposes. Its unique capability to load two carbon fibers, each with a diameter of 0.8 mm, enables the deposition of thicker carbon films when necessary or as required for specific applications.

Desk Sputter Coater

Touch Screen Control with Colorful Display

The DSCR, also known as the Desk SEM Coater, is outfitted with a 7-inch color touchscreen that offers complete automatic control and data input capabilities. This user-friendly interface is designed to be operable even by individuals with limited experience. Users can conveniently monitor vacuum levels, current readings, and deposition information, which can be displayed either as digital data or in graphical curve formats directly on the touchscreen. Additionally, the system has the capacity to store information related to the last 300 coatings, providing easy access through the history page for reference.

Desk Sputter Coater

Sample Stages

The DSCR offers versatility in sample stage configurations to cater to varying user requirements. These sample stages are designed to be rotatable, with adjustable height settings, and they can be easily interchanged. Notably, the rotary planetary sample stage is an excellent choice for achieving uniform coating, particularly when dealing with porous specimens.

Desk Sputter Coater

Desk Sputter Coater

Desk Sputter Coater

Desk Sputter Coater

THEORY & METHOD

The Desk Sputter and Carbon Coater is equipped with a rotary pump and can be configured for DC sputtering and carbon fiber coating (thread) with interchangeable heads, all within one instrument. The DSCR is an adaptable machine designed for preparing samples for scanning electron microscopy (SEM).

To enable the observation of non-conductive or poorly conductive specimens using a scanning electron microscope, their surface conductivity needs to be increased during the sample preparation phase before SEM microscopy.

Both sputtering deposition and carbon evaporation are standard methods for sample preparation in electron microscopy. The full automation option makes the system easy to use, even for operators with minimal experience. The user-friendly software with graphical interfaces simplifies the understanding of deposition information, enhancing the overall usability of the system.


TECHNICAL SPECIFICATIONS

- Two-stage, direct-drive rotary vane pump with a pumping capacity of 4m³/h
- Vacuum chamber made of Pyrex with dimensions of 170 mm outer diameter (OD) x 140 mm
- Overall dimensions: 45 cm (height) x 50 cm (width) x 37 cm (depth)
- Equipped with a 2-inch magnetron cathode
- Ultimate Vacuum: Less than 50 millitorr
- Sample stage with planetary rotation capability
- Includes a thermocouple vacuum gauge
- Quartz crystal thickness monitor with a precision of 1 nanometer
- Electrical control shutter for precise control of coating process
- Automatic venting valve for operational convenience
- Capability to save data from the last 300 coatings carried out
- Power requirements: 220V-240V, 50/60Hz, 10A

- 80 Watt DC switching power supply
- Data transfer capability via USB port to PC
- Shipping Weight: 42 kilograms

Options and Accessories:

- Carbon rod evaporation head
- High current power supply designed for carbon rod
- Quartz crystal sensor
- Planetary sample rotation capability
- Spare glass vacuum chamber
- Sputtering targets
- Carbon fiber for coating
- Sealing gaskets

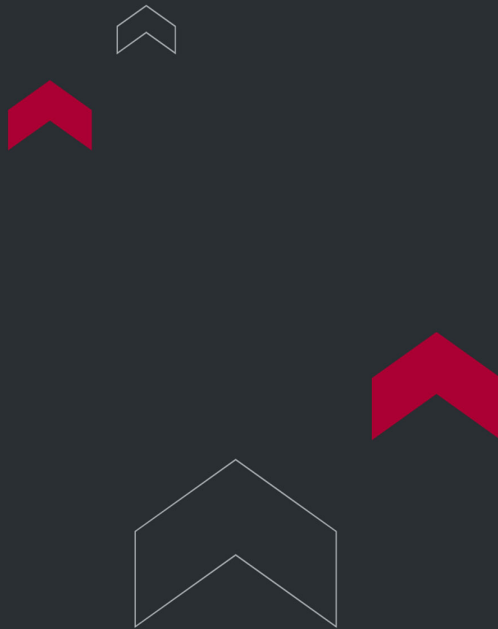

Torontech Inc.

251 Consumers Road, Suite 1200
Toronto, Ontario, M2J 4R3, Canada 

Phone: +1 416 368 2721 | Fax: +1 416 981 7652
Email: info@torontech.com

Torontech, LLC

601 Brickell Key Drive, Suite 700
Miami, FL 33131, USA 

Toll-Free: 1-866-383-7919
Email: sales@torontech.com

Torontech ME FZE

Meydan Grandstand, 6th Floor,
Meydan Road, Nad Al Sheba,
Dubai-261440, UAE 

Phone: +971 488 19252
Email: me@torontech.com

